


BSMS SHIM keypad in TopSpin

Av-300 & Av-600

BSMS Keyboard KBC

Figure 15: Layout of the BSMS Keyboard (Version BOSS)

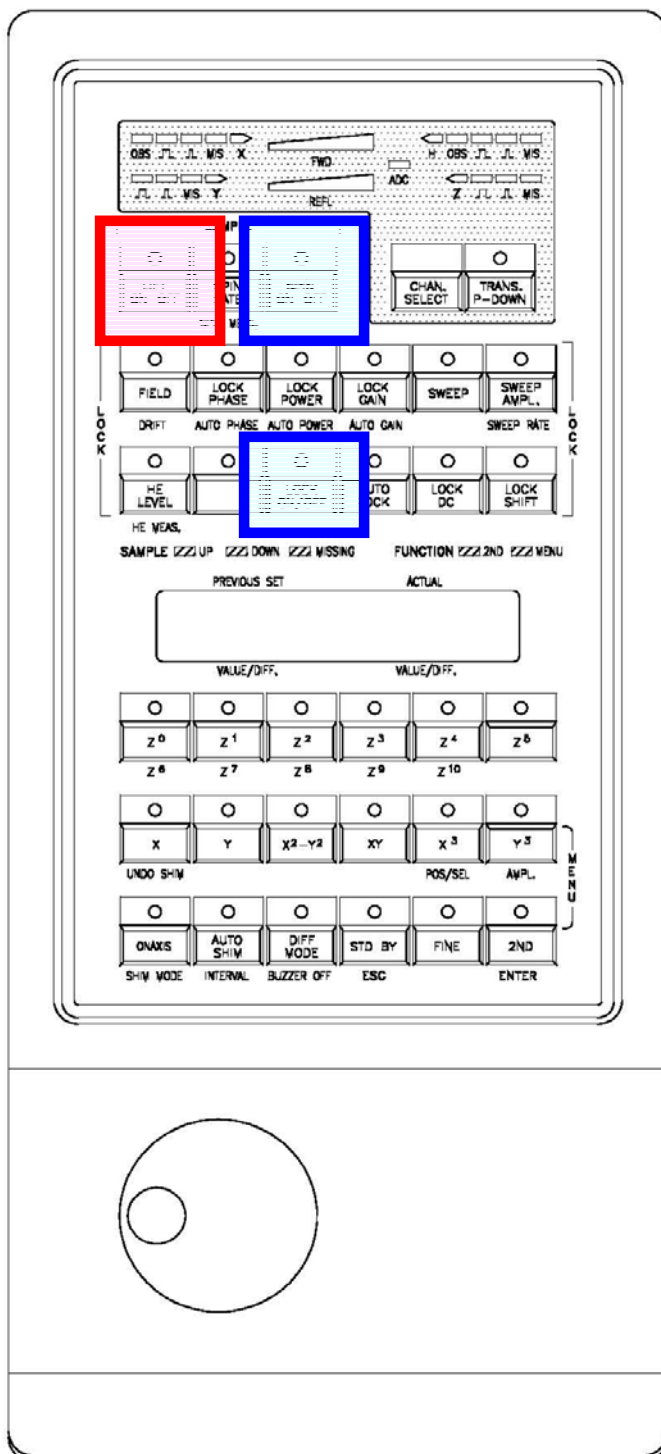

▶ halt

{if data acquisition, is in progress}

1. LOCK 'OFF'
2. SPIN 'OFF'

3. Remove bore cap

4. LIFT 'ON'

5. Change samples

6. LIFT 'OFF'

Use steps if needed when changing samples.

The magnet dewar has anti-vibration tires/legs. Do not lean/push on the side of the magnet.

Use Clean high quality NMR tubes.

Sample volume required = 0.5ml