


BSMS SHIM keypad in TopSpin

DRX-400 & DPX-300

Figure 16: Layout of the BSMS Keyboard (Version HR-20)

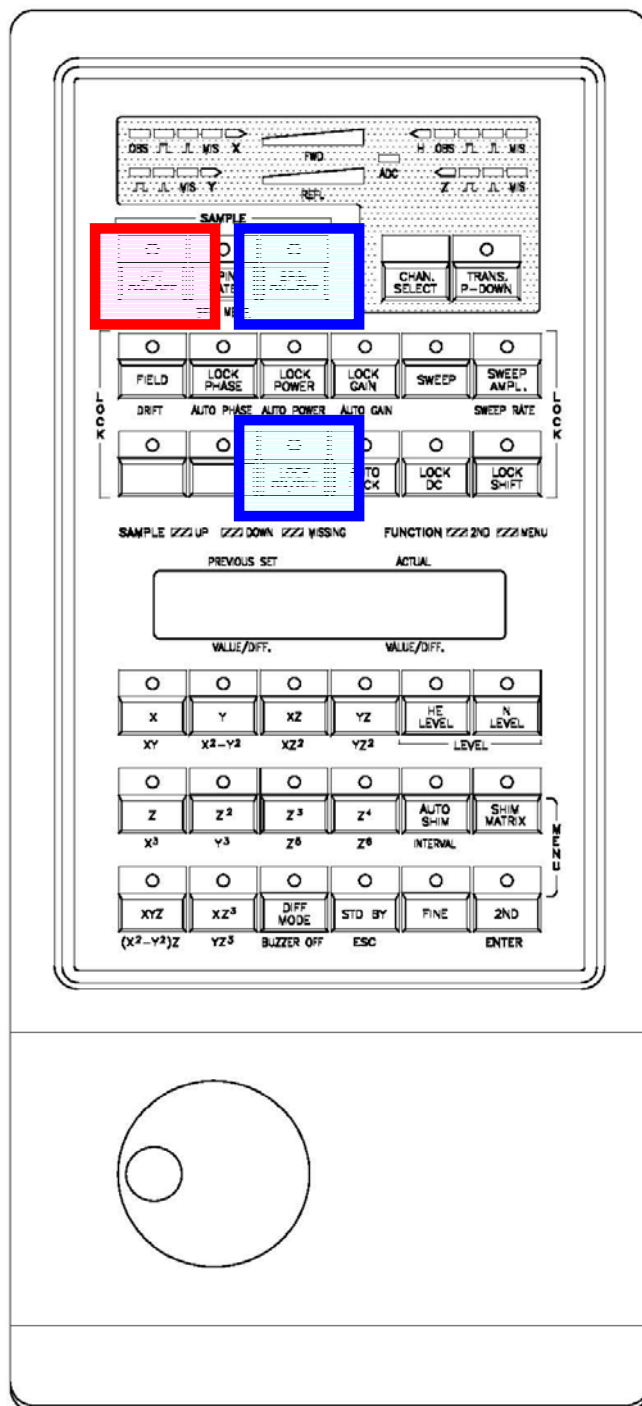

▶ **halt**
{if data acquisition, is in progress}

1. **LOCK** 'OFF'
2. **SPIN** 'OFF'

3. *Remove bore cap*

4. **LIFT** 'ON'
5. *Change samples*

6. **LIFT** 'OFF'

Use steps if needed when changing samples.

The magnet dewar has anti-vibration tires/legs. Do not lean/push on the side of the magnet.

Use Clean high quality NMR tubes.

Sample volume required = 0.5ml