

Finish . your NMR Session

TopSpin




re 1H

Click on the 're 1H' button

re 1H




Close any extra open data windows.
.....Keep **1H** data window open.

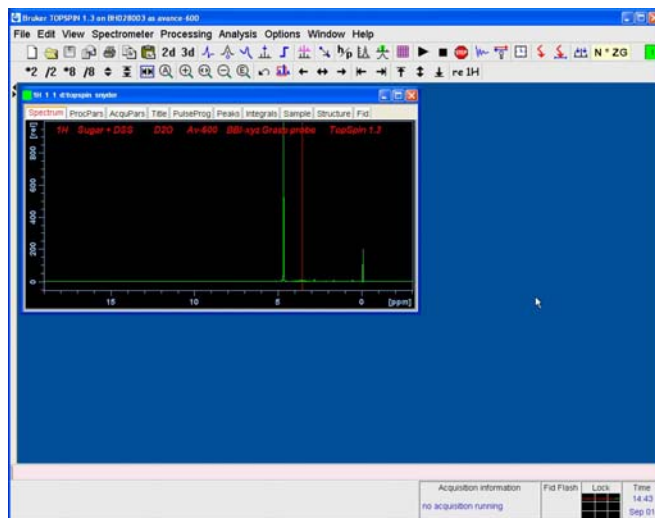



.....Ready for next TopSpin user.....

Finish . Exit TopSpinTo go to XwinNMR or to Logout.....

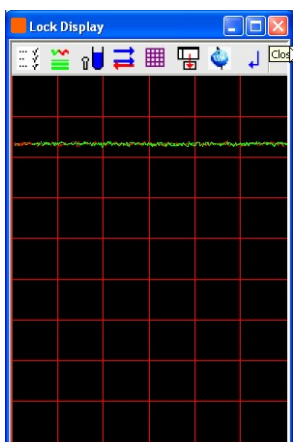

Close **Lock Display** window.

[Bruker TopSpin 1.3]
File.....
.....Exit

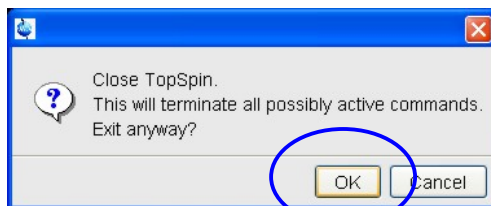
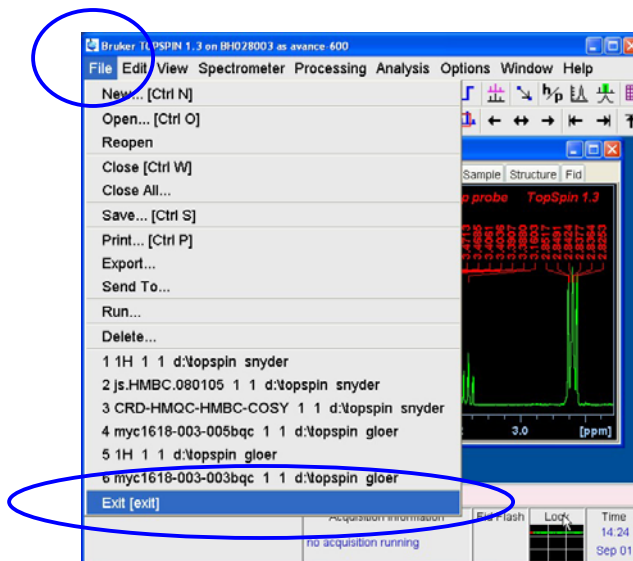


.....Ready for XwinNMR_{3.5} or TopSpin_{1.3} usage.....


Finish Logout.....



Exit from all running programs

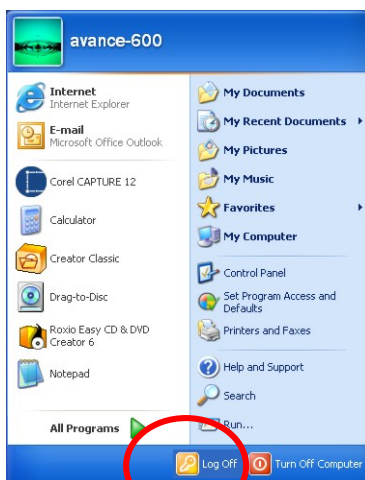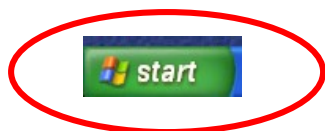

Log Off

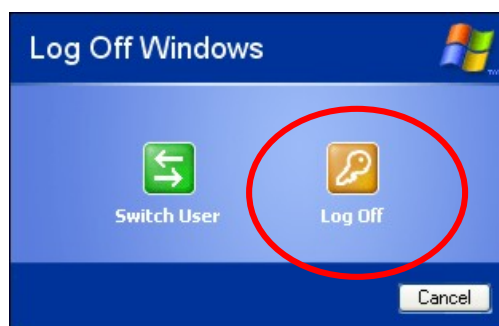

Log Off