



Wobb to Tune ^{13}C then ^1H in TopSpin

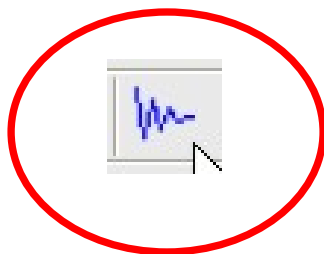

Open Acquisition Display Window [acqu]

wobb

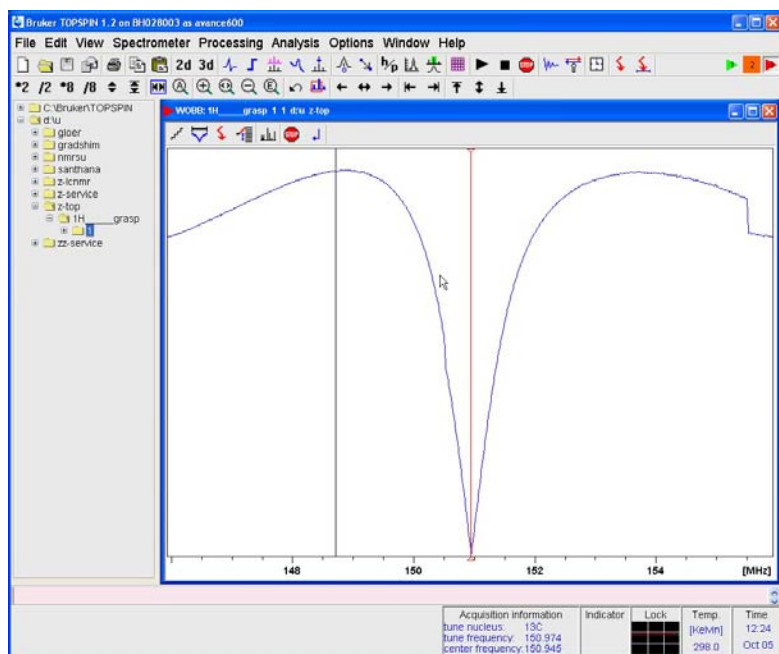


Adjust ^{13}C Tuning & Matching if needed.

Frequently for the 600 BBI-grasp probe, the DIP will be centered & no further changes are needed.



Click to switch to ^1H tuning

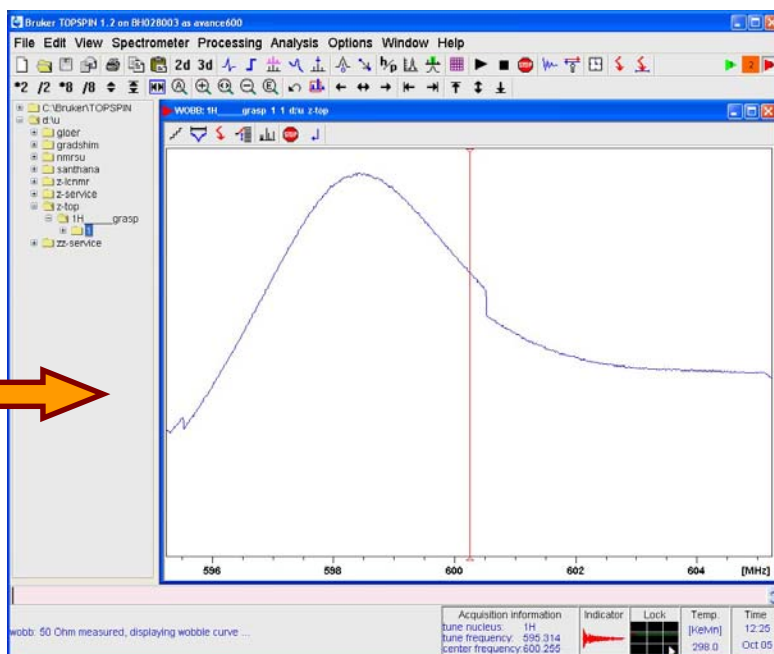


Wobb continued....

Mouse-1, and drag TopSpin Display window
to the right,

such that it is now displayed on the **2nd monitor**
which is in the **magnet-pit**.

Adjust ¹H Tuning & Matching

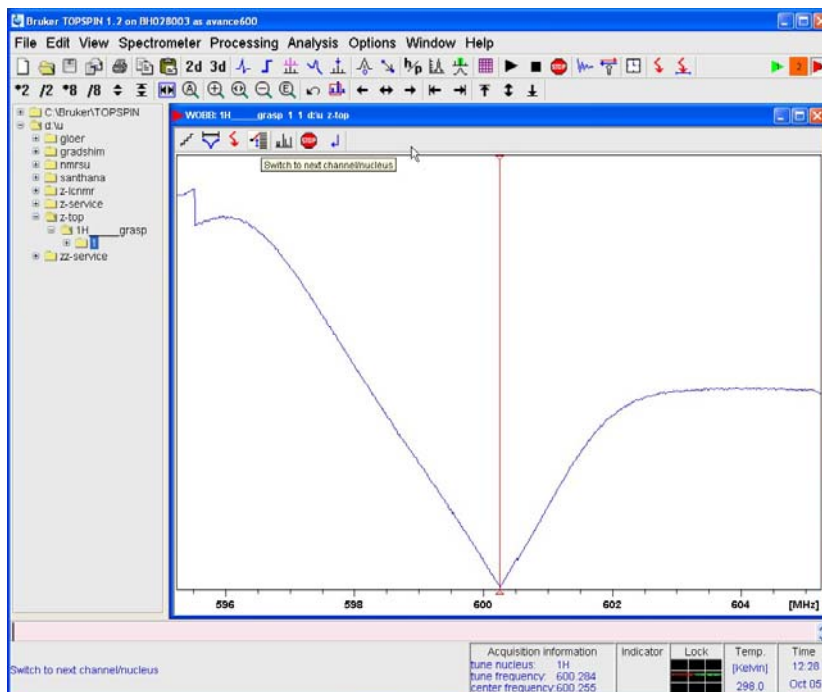

stop

Av-600

Error message that may
appear, but is of no concern.

Can ignore this message.

